
CELESTRON
Kids
Microscope
 with Smartphone Adapter
 #44119

**QUICK
 SET-UP
 GUIDE**



ENGLISH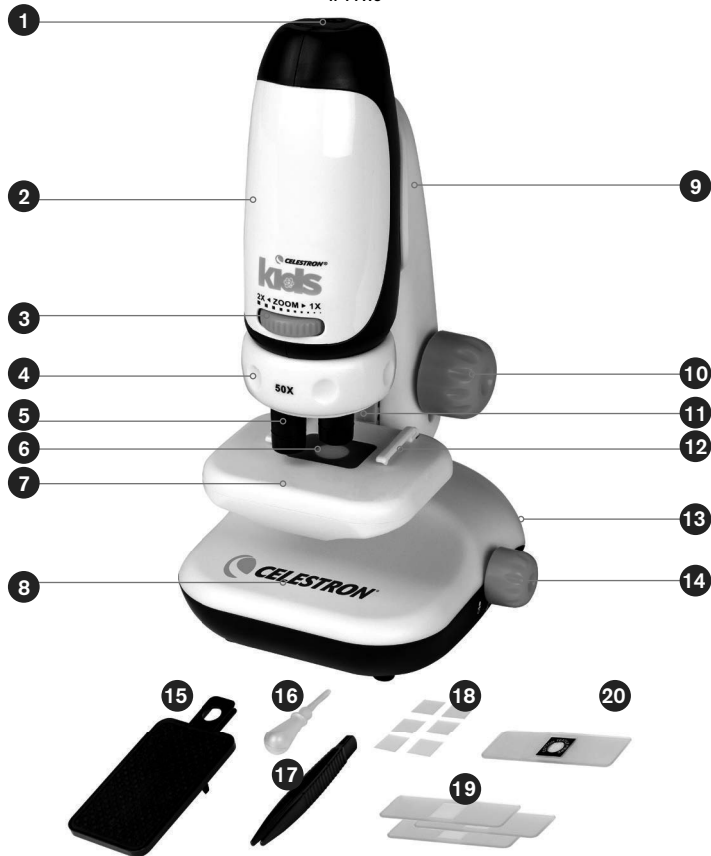

- | | |
|----------------------------|---------------------------------------|
| 1. Eyepiece / Adapter Slot | 11. Reflective LED Light |
| 2. Handheld Microscope | 12. Specimen Slide Clips |
| 3. Zoom Dial | 13. Mini Storage Drawer |
| 4. Turret | 14. LED On/Off Switch |
| 5. Objective Lens | 15. Smartphone Adapter |
| 6. Transmissive LED Cover | 16. Dropper |
| 7. Specimen Stage | 17. Tweezers |
| 8. Microscope Base | 18. 6 Blank specimen slide covers |
| 9. Microscope Arm | 19. 3 Blank specimen slides |
| 10. Focusing Dial | 20. Prepared slide with cotton swatch |



Installing the Battery

1. Open the Battery Compartment Door at the bottom of the Microscope Base (8) by lifting it up while pressing the Compartment Door Latch in the direction shown in the following diagram.
2. Install 3 fresh AA batteries with their positive (+) and negative (-) poles correctly positioned.
3. Replace the Battery Compartment Door. The microscope is now ready for use. Be sure to remove the batteries before storing the Microscope for any length of time.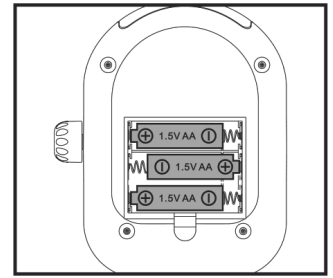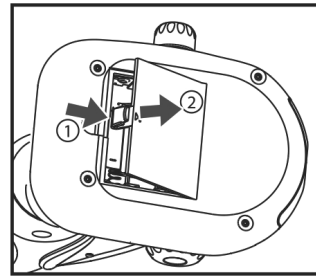




Using the Microscope

1. Place the Specimen Slide on the Specimen Stage (7) and secure with the Specimen Slide Clips (12).
2. Turn the Turret (4) to the desired magnification. (We recommend starting with 50x and increasing the magnification as needed.)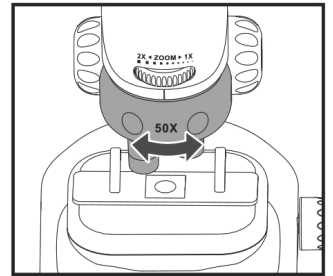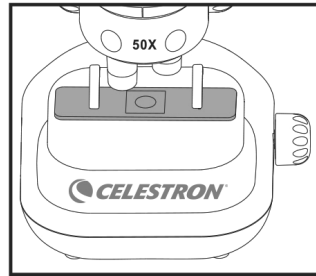

Illumination



1. Turn the LED ON/OFF Switch (14) to the right to turn on the Transmissive LED to illuminate transparent specimens.
2. Turn the LED ON/OFF Switch to the middle to turn off all LEDs
3. Turn the LED ON/OFF Switch to the left to turn on the Reflective LED (11) to illuminate non-transparent specimens.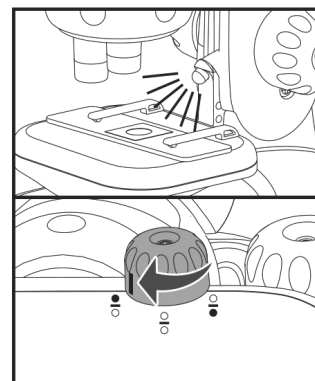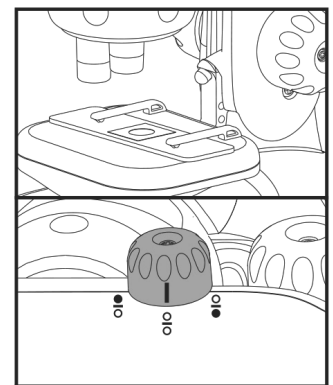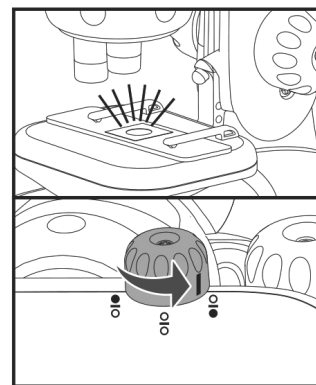

Focusing

Turn the Focusing Dial (10) clockwise to rise the Specimen Stage (7) to the highest position toward the Objective Lens (5). Take caution to ensure that the Specimen Slide does not touch the Objective Lens.

View through the eyepiece and lower the Specimen Stage gradually by turning the Focusing Dial counter-clockwise until the image is sharp and clear. You may need to move the Specimen Slide to locate the desired area of specimen examination.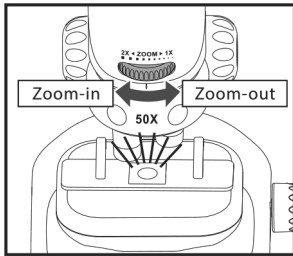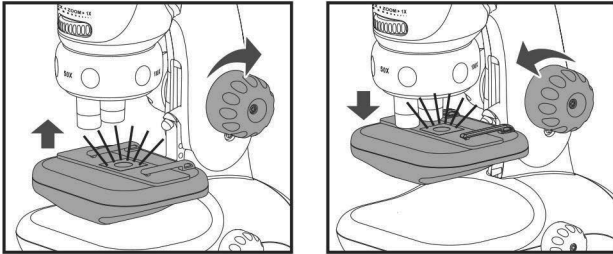


Zoom Dial

The Zoom Dial (3) may be used to zoom in on the image.

Zoom Magnification: 50x-100x, 100x-200x, 200-400x


Microscope Arm

Pull back on the Microscope Arm (9) to adjust viewing angle.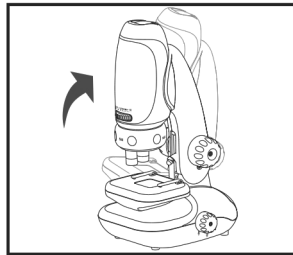

Handheld


Remove the microscope from the Microscope Arm for handheld use. Sufficient lighting is required to view image with Handheld Microscope (Not recommended for indoor use.) Always start at 50x magnification before zooming in or turning to a higher magnification.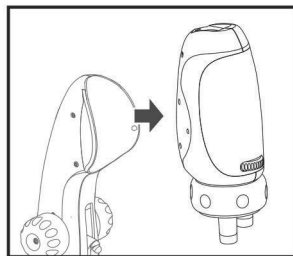

Smartphone Adapter

CAUTION: Before placing a smartphone on the smartphone adapter, check to ensure that Microscope Arm is in the upright position and that the Handheld Microscope is securely attached to the Microscope Arm to prevent it from slipping and the smartphone being damaged. Do not use the Smartphone Adapter with the Microscope Arm in any position other than upright.

1. Remove the Adapter Slot (1) cover from the microscope.
2. Slide the Smartphone Adapter (15) into the Adapter Slot (1).
3. Align the camera of the smartphone to the Eyepiece (1) and rest the smartphone on the Smartphone Adapter (15).

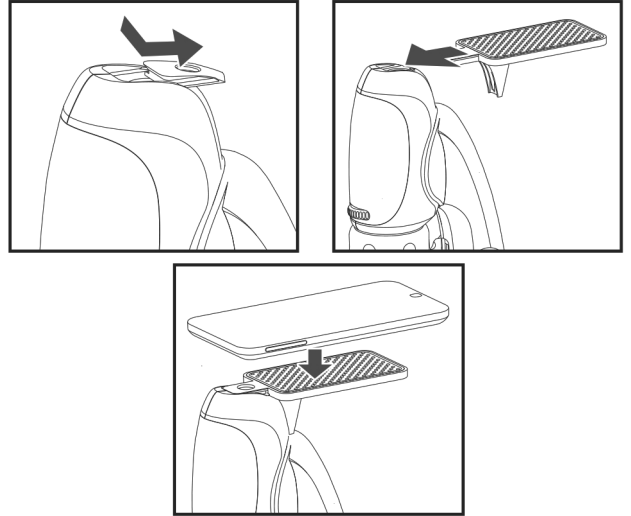Specifications

Magnifications	50x-100x, 100x-200x, 200x-400x
Optical	3-Objective Turret, 2x optical zoom
Effective Observation area (without Zoom)	2.3mm (50x), 1.2mm (100x), 0.56mm (200x)
Effective Observation area (with max zoom)	1.2mm (100x), 0.6mm (200x), 0.28mm (400x)
Power	3 x 1.5V AA alkaline batteries (not included)
Illumination	Two LEDs
Size	112mm x 137mm x 224mm
Weight	Approx. 335 grams (excluding batteries)

Safety Information

1. To avoid possible permanent eye damage, do not look at the Sun or bright objects with the microscope.
2. To avoid possible permanent eye damage, do not place the lighted microscope directly upon the human eye.
3. Remove the battery before storing the microscope for an extended period.
4. Do not short-circuit the supply terminals.
5. Do not mix different types of batteries or new and used batteries.
6. Do not recharge nonchargeable batteries.
7. Remove rechargeable batteries from the device before being charged.
8. Only charge rechargeable batteries under adult supervision.
9. Remove exhausted batteries from the device.
10. Please retain the packaging and instruction manual for future reference.

CHOKING HAZARD! This kit contains small parts. Not suitable for children under 3 years.

WARNING: Never point your Microscope at the Sun. It will cause serious damage. Always use your Microscope under adult supervision.

© 2020 Celestron • All rights reserved • celestron.com/pages/technical-support
2835 Columbia Street, Torrance, CA 90503 U.S.A. • Made in China • This product is designed and intended for use by those 8 years of age and older.

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