ENGLISH

8. **ROTATE THE FOCUS WHEEL [B]** to bring object in clear focus.
9. **PRESS THE CAPTURE BUTTON** to take a photo or record video.
10. **ZOOM IN ON OBJECT** using the up arrow. Zoom out using the down arrow.

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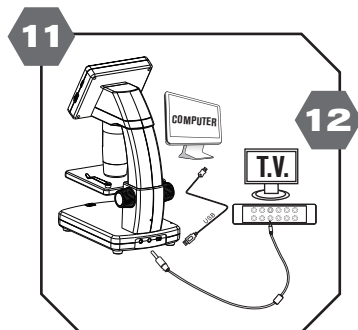
11. **CONNECT THE MICROSCOPE** to the computer by the included USB cable to A.) Transfer Files or B.) Preview image on computer screen.
12. **CONNECT THE MICROSCOPE** to a TV monitor by A.) inserting the 1/8" plug on the end of the supplied AV cable into the microscope Video port (on back of base) and the RCA plug into the yellow "video" input connector on TV.

NOTE: For full manual instructions go to the product page at celestron.com

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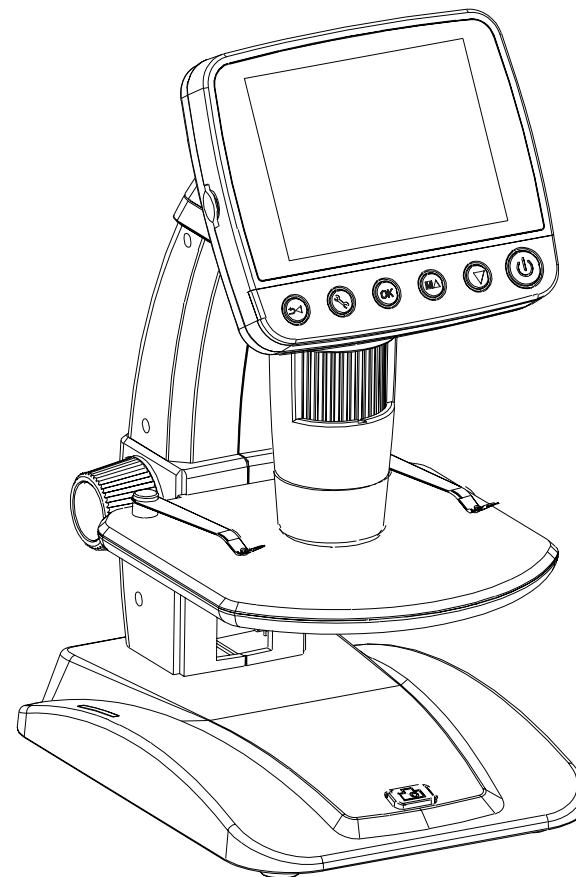
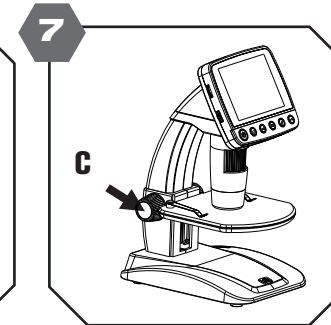
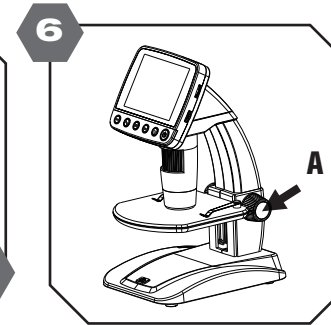
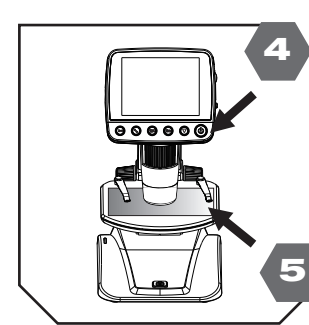
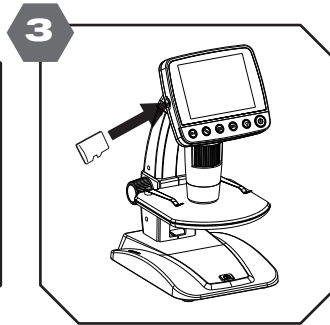
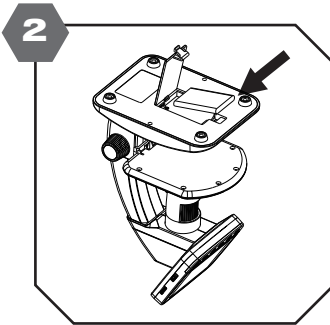
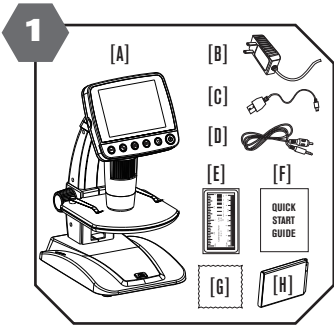
2. INSTALL BATTERY

- [1] Open Battery door.
- [2] Insert Lithium Battery following the polarities shown inside the compartment.

- 3. **INSERT MICRO SD CARD**** into the micro SD card slot; gently press in until card is latched.

NOTE: You can still view images on LCD screen without the Micro SD card, they cannot be saved to the unit. See manual on CD for complete instructions on using the software and your PC.

- 4. **TURN ON MICROSCOPE** by pressing On/Off button.
- 5. **PLACE OBJECT** to be observed on the stage.
- 6. **ROTATE THE STAGE ADJUSTMENT WHEEL [A]** to bring the object near the microscope tube.
- 7. **LOCKING THE STAGE IN PLACE [C]** by rotating the stage locking wheel.

*44361 - AC Multi-Country Plug Adapter (1) **Micro SD Card Not Included ***
NOTE: You can still view images on LCD screen without the Micro SD card, they cannot be saved to the unit.

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